

AC600 Semi-Automated Calorimeter

Specification Sheet



| | | |
|--|--|--|
| Method | Isoperibol | |
| Range | | |
| Joules/Charge: | 14,000 [†] to 35,000 ^{††} (14,000 [†] to 35,000 ^{††} kJ/kg for a 1 gram sample) | |
| Calories/Charge: | 3,300 [†] to 8,300 ^{††} (3,300 [†] to 8,300 ^{††} kcal/kg for a 1 gram sample) | |
| BTU/Charge: | 13 [†] to 33 ^{††} (6,000 [†] to 15,000 ^{††} BTU/lb for a 1 gram sample) | |
| Precision | 0.10% RSD* | |
| Analysis Time | TruSpeed [®] Mode: 5 minutes | Delta T Mode: 9 minutes |
| Corrections | Nitrogen, sulfur, fuse wire, moisture, spike, and ash | |
| User-Selectable Ignition | String or wire-fuse | |
| Temperature Measuring Resolution | 0.0001°C | |
| Gas Requirements | Analytical: Oxygen 450 psi (31.0 bar) max; 99.5% purity Pneumatic: Compressed Air; 12 psi (0.8 bar); <i>source must be oil and water free</i> | |
| Electrical Requirements | 90 to 127 V~/180 to 260 V~(at max load), 50/60 Hz, single phase, 5 A, 2,000 BTU/hr | |
| Water Requirements | | |
| Recirculating Water Chiller [‡] Nominal Temp: | 15°C (59°F) | |
| Cooling Capacity (at 15°C): | Single AC600: 680 BTU/hr | Dual AC600: 1,400 BTU/hr |
| Type: | Distilled Water <i>NOTE: Do NOT use deionized water</i> | |
| pH: | 6-8 | |
| Dissolved Solids: | 0.5 to 100 ppm | |
| Resistivity: | 50 kΩ•cm to 2 MΩ•cm (0.5 μs/cm to 20 μs/cm) | |
| Pressure: | 4 to 15 psi (0.3 to 1.0 bars) | |
| Volume: | Single AC600: 3 gallons (12 liters) for initial setup Dual AC600: 9 gallons (35 liters) for initial setup | |
| Required Work Space** | 15 in. W x 21 in. D x 18 in. H (38 cm x 53 cm x 46 cm) | |
| Weight (approx.) | 135 lb. (61 kg) | Shipping Weight (approx.): 148 lb. (67 kg) |
| Environmental Conditions | Operating Temp: 15° to 35°C (59° to 95°F) | Humidity: 20 to 80%, non-condensing |
| | General: Area should be free of drafts, and shielded from sunlight and other radiation sources | |
| Part Numbers | | |
| AC600SC | AC600 Semi-Automatic Calorimeter with standard vessel; vessel preparation station; external PC | |
| AC600SHC | AC600 Semi-Automatic Calorimeter with halogen-resistant vessel; vessel preparation station; external PC | |
| AC600DC | AC600 Dual Semi-Automatic Calorimeter with two standard vessels; vessel preparation station; external PC | |
| AC600DHC | AC600 Dual Semi-Automatic Calorimeter with two halogen-resistant vessels; vessel preparation station; external PC | |

*Based on analysis of benzoic acid at 1 gram; n = 7.

**Allow a 6-inch (15 cm) minimum access area around instrument.

†Lower values can be measured by spiking samples that are not completely combusting.

††This is the combustion vessel safety limit. Do NOT exceed this limit. Exceeding this limit could result in vessel failure causing death, serious personal injury, and/or property damage.

‡Chiller must be capable of operating continuously with flow stopped to the AC600.

V~ denotes VAC.


Theory of Operation

The LECO AC600 semi-automatic calorimeter is a Windows®-based instrument developed to measure the calorific content of various organic materials such as coal, coke, and fuel oil.

The calorific value of a sample is determined by precisely measuring the heat released after combustion of the sample in a controlled environment. The heat released is proportional to the calorific value of the substance.

The sample is placed into a combustion vessel, which is pressurized with oxygen. The combustion vessel is automatically lowered into a water bath within the instrument known as a bucket. The bucket is sealed using an inflatable bladder. The cavity surrounding the bucket is known as the jacket and is also filled with water. The water temperature in the jacket is closely controlled at a precise set temperature (isoperibol calorimeter system). The sample is ignited and the temperature of the bucket and jacket water is measured by an electrical thermometer with a resolution of 0.0001 of a degree. A measurement of the water temperature inside the bucket and jacket is collected every second. The results can be corrected within the software for any spiking, nitrogen, sulfur, moisture, and ash content if necessary. Two options for analyzing data are available. A simple temperature difference may be chosen by the user (Delta T mode), in which calorific values are determined by a simple maximum temperature rise of the bucket. The user may also choose to use a thermodynamic model (TruSpeed mode) that has been developed by LECO to model heat exchange within the AC600 system. The TruSpeed mode takes into account heat capacities of the system's components along with corrections for energy transfer within the system. The TruSpeed mode enables the AC600 to attain rapid 5 minute analysis times without compromising the accuracy or precision of the calorific result.

Flow Diagram

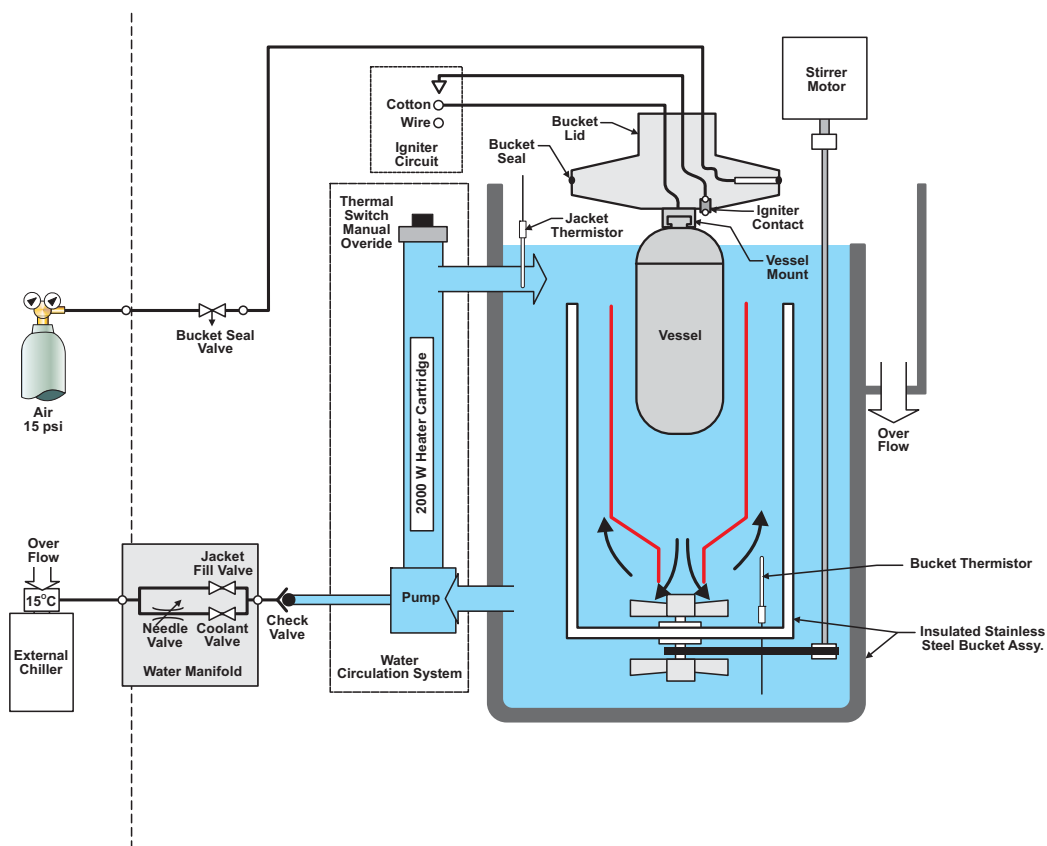

Specifications and part numbers may change.
Consult LECO for latest information.

LECO, TruSpeed are registered trademarks of LECO Corporation.
Microsoft, Windows are registered trademarks of Microsoft Corporation.

3000 Lakeview Avenue • St. Joseph, MI 49085 • Phone: 800-292-6141 • Fax: 269-982-8977
info@leco.com • www.leco.com • ISO-9001:2008 HQ-Q-994

LECO Corporation



Delivering the Right Results

© 2015 LECO Corporation